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### 1. Introduction

These general instructions must be readily available to operators of the electrode at all times. Before using the electrode, these instructions must be read and understood by all operating personnel.

The InPro 4501 VP electrode is supplied with a VarioPin connector. The VarioPin connector is used to connect the electrode to the transmitter.

The METTLER TOLEDO electrode versions InPro 4501 VP and InPro 4501 fixed cable are low-maintenance, temperature-compensated pH combination electrodes containing polymer electrolyte (XEROLYT® PLUS), and are hereafter referred to as the InPro 4501 family. An integrated temperature sensor (RTD) provides automatic compensation of the electrode slope to ensure accurate readings. The InPro 4501 family is suitable for pH measurements in:

- emulsions and suspensions
- sulfide-bearing media
- solutions containing heavy metals (lead, mercury, silver, etc.)
- processes with rapid pressure fluctuations (within the given range)
- measurements in wastewater and in decontamination processes for industrial effluents
- heavily contaminated media

The InPro 4501 family electrodes are unsuitable for sterilization/autoclaving procedures, for processes with rapid temperature fluctuations or food and beverage applications.

The InPro 4501 VP electrode is supplied with a VarioPin connector. The VarioPin connector is used to connect the electrode to the transmitter.

### 2. Safety remarks

The InPro 4501 family of electrodes is intended solely for the purpose described in §1 above. The electrodes consist partly of glass. If they are to be cleaned or calibrated using acid or alkaline solutions, protective clothing (goggles and gloves) must be worn. In addition, local and in-house-safety regulations must be observed. We recommend that the electrodes only be put into operation in conjunction with original products/accessories from METTLER TOLEDO. The electrodes may be operated and serviced only by personnel familiar with the equipment in question and who have read and understood these instructions.

The InPro 4501 VP electrode is supplied with a VarioPin connector. The VarioPin connector is used to connect the electrode to the transmitter.

### 3. Product description

The markings on each electrode provide the following information:

InPro®4501VP/xxx/yyy	designation / temperature sensor / solution ground (SG)
InPro®4501/xxx/yyyy/zzz/www	fixed cable designation/temperature sensor/solution ground (SG)/cable length /cable end
combination electrode	type of sensor
Pt100 / Pt1000	type of built-in temperature sensor
pH 1...14 / 0...100 °C	pH measuring range / temperature range
Order no. 41 453 310x	ordering number VP
Order no. 10 453 30xx	ordering number fixed cable

### 4. Installation and preparation for use

- On unpacking, and again prior to installation, check the electrode visually for signs of mechanical damage to the glass membrane, the shaft and the connector. Report any faults or defects immediately to your authorized METTLER TOLEDO dealer.
- Remove the watering cap and rinse the electrode with distilled or de-ionized water. Use only soft tissue to dab the electrode dry. Do not rub the glass membrane since this can lead to sluggish response behavior of the electrode.
- Electrodes with a dry glass membrane must be conditioned for several hours in a 3mol KCl solution before being used.

- Connect always the solution ground (blue cable) to the transmitter.
- Use Teflon tape to seal the NPT thread.
- If possible use the protection sleeve (order no. 52 401 808).

## 5. Cabling

The InPro 4501 VP is equipped with a VarioPin connector. To connect the cable to the electrode, attach and gently rotate until the connector is engaged (groove). Screw the cable cap nut onto the electrode until the connection is secure. The plug assembly should be fitted hand-tight only; no tools should be used. For details of the exact assignment of the individual cable leads to the VP pins, please refer to the instruction leaflet "Preparation of the cable for connection to the transmitter" supplied together with the cable.

The InPro 4501 VP electrode is supplied with a VarioPin connector. The VarioPin connector is used to connect the electrode to the transmitter.

InPro4501 fixed cable. There are two options for termination for the InPro 4501 with fixed cable:

- BNC connector for pH combination electrode with stripped and tinned leads for RTD and solution ground.
- Stripped and tinned leads for pH combination electrode, for RTD and solution ground.

Refer to your transmitter instruction manual for correct wiring.

**NOTE:** The solution ground lead is designed to go to a galvanically isolated terminal on the transmitter and should not be used as an outer ground shield. If the transmitter does not have sensor diagnostics, this lead should be left unterminated.

<b>Color</b>	<b>Function</b>
black/transparent	glass
red	coax shield/reference
grey	RTD sensing (temperature sensor, 3-wire connection)
blue	solution ground
white	RTD return (temperature sensor, 2-wire connection)
green	RTD in (temperature sensor, 2-wire connection)
green/yellow	cable shield

## 6. Calibration and measurement

InPro4501VP electrodes require two-point calibration. Before calibration, the watering cap must be removed from the pH electrode. The electrode is then dipped successively into two different buffer solutions, each of known pH value, and the transmitter calibrated to these buffer values (for example, buffers pH 4.01 and pH 7.00). Please note the instructions given in the operating manual of the transmitter. Calibration should be repeated as and when needed.

The InPro 4501 VP electrode is supplied with a VarioPin connector. The VarioPin connector is used to connect the electrode to the transmitter.

### 7. Maintenance

- Electrodes with XEROLYT®PLUS polymer electrolyte cannot be refilled.
- Electrodes should not be stored dry for any period of time
- During storage, the watering cap should be in place, filled with FRISCOLYT or 3mol KCl solution.
- Never store an electrode in distilled or deionized water, as this shortens the lifetime of the electrode.
- Should the electrode dry out, it must be conditioned for 24 hours in FRISCOLYT (ordering no. 51 2340 053 for 250 ml, or 51 340 054 for 6 x 250 ml) before reuse. After conditioning, recalibration is required.
- After each workcycle, the tip of the electrode should be cleaned using de-ionized water.

**ATTENTION:** Do not under any circumstance use METTLER TOLEDO reactivating solution (ordering no. 51 319 053 / in USA 43 989 5125) as this contains hydrofluoric acid (HF) which would attack the XEROLYT reference system.

The InPro 4501 VP electrode is supplied with a VarioPin connector. The VarioPin connector is used to connect the electrode to the transmitter.

### 8. Diagnostics

The electrodes are equipped with a grounding ring which allows a diagnostic of the diaphragm condition. If this diaphragm (open junction) should become clogged or if the reference resistance exceeds a certain value (see transmitter operation manual), it is then necessary to replace the electrode. Electrodes of the InPro 4501 family should as a general rule not be cleaned with acidic cleaning solutions, e.g. diaphragm cleaning solution, reactivation solution, etc. However, by extremely careful application, it is possible to reactivate the glass membrane, see «point 9» below.

The InPro 4501 VP electrode is supplied with a VarioPin connector. The VarioPin connector is used to connect the electrode to the transmitter.

### 9. Troubleshooting (sluggish response time, insufficient sensitivity, etc.)

**ATTENTION:** The glass membrane must be very cautiously dabbed with HF solution using a fine brush and taking care that solution does not come into contact with the titanium pin (metallic solution ground). Also check the electrode connector (only VP version) for possible traces of moisture. If necessary, clean well using deionized water or alcohol, and afterwards dry carefully.

The InPro 4501 VP electrode is supplied with a VarioPin connector. The VarioPin connector is used to connect the electrode to the transmitter.

### 10. Disposal

Dispose of InPro 4501 family electrodes in accordance with local or state regulations.

The InPro 4501 VP electrode is supplied with a VarioPin connector. The VarioPin connector is used to connect the electrode to the transmitter.

### 11. Guarantee

METTLER TOLEDO grants 12 months warranty covering manufacturing defects, valid from the date of production of the electrode.