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General instructions for pH combination electrode InPro®4550VP

1. Introduction

These general instructions must be readily available to operators of the electrode at all times. Before using the electrode, these instructions must be read and understood by all operating personnel.

The METTLER TOLEDO electrode versions InPro 4550VP are low-maintenance, temperature-compensated pH combination electrodes containing polymer electrolyte (XEROLYT® PLUS), and are hereafter referred to as the InPro 4550 family. An integrated temperature sensor (RTD) provides automatic compensation of the electrode slope to ensure accurate readings.

The InPro 4550 family is suitable for pH measurements in:

- emulsions and suspensions
- sulfide-bearing media
- solutions containing heavy metals (lead, mercury, silver, etc.)
- processes with rapid pressure fluctuations (within the given range)
- heavily contaminated media.

The InPro 4550 family electrodes are unsuitable for sterilization/autoclaving procedures, for processes with rapid temperature fluctuations or food and beverage applications.

2. Safety remarks

The InPro 4550 family of electrodes is intended solely for the purpose described in §1 above. The electrodes consist partly of glass. If they are to be cleaned or calibrated using acid or alkaline solutions, protective clothing (goggles and gloves) must be worn. In addition, local and in-house-safety regulations must be observed. We recommend that the electrodes only be put into operation in conjunction with original products/accessories from METTLER TOLEDO. The electrodes may be operated and serviced only by personnel familiar with the equipment in question and who have read and understood these instructions.

3. Product description

The markings on each electrode provide the following information:

InPro®4550VP/xxx/yyy	designation / temperature sensor / solution ground (SG)
combination electrode	type of sensor
Pt100 / Pt1000	type of built-in temperature sensor
pH 0...14 / 0...130 °C	pH measuring range / temperature range
Ⓢ II 1/2G EEx ia IIC T6/T5/T4	Ex-classification
SNCH 00 ATEX 3130X	designation and number of certificate
IS CL I, Div1, GR ABCD	FM approved
Order no. 52 002 40x	ordering number

4. Installation and preparation for use

- On unpacking, and again prior to installation, check the electrode visually for signs of mechanical damage to the glass membrane, the shaft and the connector. Report any faults or defects immediately to your authorized METTLER TOLEDO dealer.
- Remove the watering cap and rinse the electrode with distilled or de-ionized water. Use only soft tissue to dab the electrode dry. Do not rub the glass membrane since this can lead to sluggish response behavior of the electrode.
- Electrodes with a dry glass membrane must be conditioned for several hours in a 3mol KCl solution before being used.
- Connect always the solution ground (blue cable) to the transmitter.
- Use Teflon tape to seal the NPT thread.
- If possible use the protection sleeve (order no. 52 401 808).

5. Cabling

The InPro 4550VP is equipped with a VarioPin connector. To connect the cable to the electrode, attach and gently rotate until the connector is engaged (groove). Screw the cable cap nut onto the electrode until the connection is secure. The plug assembly should be fitted hand-tight only; no tools should be used. For details of the exact assignment of the individual cable leads to the VP pins, please refer to the instruction leaflet «Preparation of the cable for connection to the transmitter» supplied together with the cable.

6. Calibration and measurement

InPro 4550VP electrodes require two-point calibration. Before calibration, the watering cap must be removed from the pH electrode. The electrode is then dipped successively into two different buffer solutions, each of known pH value, and the transmitter calibrated to these buffer values (for example, buffers pH 4.01 and pH 7.00). Please note the instructions given in the operating manual of the transmitter. Calibration should be repeated as and when needed.

7. Maintenance

- Electrodes with XEROLYT®PLUS polymer electrolyte cannot be refilled.
- Electrodes should not be stored dry for any period of time.
- During storage, the watering cap should be in place, filled with FRISCOLYT or 3mol KCl solution.
- Never store an electrode in distilled or deionized water, as this shortens the lifetime of the electrode.
- Should the electrode dry out, it must be conditioned for 24 hours in FRISCOLYT (ordering no. 51 2340 053 for 250 ml, or 51 340 054 for 6 x 250 ml) before reuse. After conditioning, recalibration is required.
- After each workcycle, the tip of the electrode should be cleaned using de-ionized water.

ATTENTION: Do not under any circumstance use METTLER TOLEDO reactivating solution (ordering no. 51 319 053 / in USA 43 989 5125) as this contains hydrofluoric acid (HF) which would attack the XEROLYT reference system.

8. Diagnostics

The electrodes are equipped with a grounding ring which allows a diagnostic of the diaphragm condition. If this diaphragm (open junction) should become clogged or if the reference resistance exceeds a certain value (see transmitter operation manual), it is then necessary to replace the electrode. Electrodes of the InPro 4550 family should as a general rule not be cleaned with acidic cleaning solutions, e.g. diaphragm cleaning solution, reactivation solution, etc. However, by extremely careful application, it is possible to reactivate the glass membrane, see point 9 below.

9. Troubleshooting (sluggish response time, insufficient sensitivity, etc.)

ATTENTION: The glass membrane must be very cautiously dabbed with HF solution using a fine brush and taking care that solution does not come into contact with the titanium pin (metallic solution ground). Also check the electrode connector (VP) for possible traces of moisture. If necessary, clean well using deionized water or alcohol, and afterwards dry carefully.

10. Disposal

Dispose of InPro 4550 family electrodes in accordance with local or state regulations.

11. Guarantee

METTLER TOLEDO grants 12 months warranty covering manufacturing defects, valid from the date of production of the electrode.

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Subject to technical changes. 01/06 © Mettler-Toledo GmbH. Printed in Switzerland